

Instructions for Use

Life Science Kits & Assays



innuPREP RNA Kit - PP Mini

Order No.:

845-PS-0060016	16 reactions
845-PS-0060096	96 reactions

Publication No.: HB_PS-0060_e_230720

This documentation describes the state at the time of publishing.
It needs not necessarily agree with future versions. Subject to change!

Print-out and further use permitted with indication of source.

© Copyright 2023, IST Innuscreen GmbH

Manufacturer and Distributor:

IST Innuscreen GmbH	Phone +49 30 9489 3380
Robert-Rössle-Straße 10	Fax +49 30 9489 3381
13125 Berlin · Germany	
Made in Germany!	info.innu@ist-ag.com

Contents

1	Introduction.....	2
1.1	Intended use.....	2
1.2	Notes on the use of this manual and the kit.....	3
2	Safety precautions.....	4
3	General notes and safety recommendations on handling RNA.....	5
4	Storage conditions	6
5	Functional testing and technical assistance.....	7
6	Product use and warranty.....	7
7	Kit components	8
7.1	Components included in the kit.....	8
7.2	Components not included in the kit	8
7.3	Components needed for isolation from bacteria (not included)	8
7.4	Components needed for isolation from yeasts (not included)	9
8	Product specifications	9
9	Initial steps before starting.....	9
10	Sample Preparation	10
10.1	Protocol 1: RNA extraction from tissue samples	10
10.2	Protocol 2: RNA extraction from eukaryotic cells	12
10.3	Protocol 3: RNA extraction from bacteria and yeasts.....	12
11	Automated extraction using PurePrep Mini	14
11.1	Prefilling of the DW Plate or the DW Strips	14
11.2	Loading filled Deep Well Plate/Strips to the PurePrep Mini and plug in the Tip Combs	14
12	Troubleshooting.....	16

1 Introduction

1.1 Intended use

The innuPREP RNA Kit – PP Mini has been designed for automated isolation of total RNA from tissue sample, eukaryotic cells, bacterial cell pellets and yeasts.

The extraction procedure takes place on the magnetic particle processor PurePrep Mini and allows the parallel and flexible extraction of up to 16 samples.

The extraction process starts with sample lysis (external step) followed by removing of DNA based on binding of DNA on a spin filter membrane. After DNA removing the filtrate is used for the automated extraction of total RNA on PurePrep Mini.

The kit is intended for use by professional users. The kit has been designed to be used for a wide range of different downstream applications, like amplification reactions and further analytical procedures.



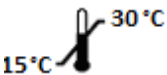







CONSULT INSTRUCTION FOR USE

This package insert must be read carefully before use. Package insert instructions must be followed accordingly. Reliability of results cannot be guaranteed if there are any deviations from the instructions in this package insert.

1.2 Notes on the use of this manual and the kit

For easy reference and orientation, the manual and labels use the following warning and information symbols as well as the shown methodology:

Symbol	Information
	REF Catalogue number.
	Content Contains sufficient reagents for <N> reactions.
	Storage conditions Store at room temperature or shown conditions respectively.
	Consult instructions for use This information must be observed to avoid improper use of the kit and the kit components.
	Expiry date
	Lot number The number of the kit charge.
	Manufactured by Contact information of manufacturer.
	For single use only Do not use components for a second time.
	Note / Attention Observe the notes marked in this way to ensure correct function of the kit and to avoid operating errors for obtaining correct results.

The following systematic approach is introduced in the manual:

- The chapters and figures are numbered consecutively.
- A cross reference is indicated with an arrow (e.g. → „Notes on the use of this manual“ p. 3).
- Working steps are numbered.

2 Safety precautions

NOTE

Read through this chapter carefully before use to guarantee your own safety and a trouble-free operation.

Follow all the safety instructions explained in the manual, as well as all messages and information, which are shown.

All due care and attention should be exercised in handling the materials and reagents contained in the kit. Always wear gloves while handling these reagents and avoid any skin contact! In case of contact, flush eyes or skin with a large amount of water immediately.



FOR SINGLE USE ONLY!

This kit is made for single use only!

ATTENTION!

Don't eat or drink components of the kit!

The kit shall only be handled by educated personnel in a laboratory environment!

If the buffer bottles are damaged or leaking, wear gloves and protective goggles when discarding the bottles in order to avoid any injuries. This kit could be used with potential infectious samples. Therefore, all liquid waste must be considered as potentially infectious and must be handled and discarded according to local safety regulation.

Please observe the federal, state and local safety and environmental regulations. Follow the usual precautions for applications using extracted nucleic acids. All materials and reagents used for DNA or RNA isolation should be free of DNases or RNases.

ATTENTION!

Do not add bleach or acidic components to the waste after sample preparation!

NOTE

Emergency medical information in English and German can be obtained 24 hours a day from:

Poison Information Center, Freiburg / Germany

Phone: +49 (0)761 19 240.

For more information on GHS classification and the Safety Data Sheet (SDS) please contact sds.innu@ist-ag.com.

3 General notes and safety recommendations on handling RNA

RNA is far less stable than DNA. It is very sensitive to degradation by endogenous RNases in the biological material and exogenous RNases which are permanently present everywhere in the lab. To achieve satisfactory qualitative and quantitative results in RNA preparations, contaminations with exogenous RNases have to be reduced to a minimum in accordance with the following recommendations:

- Always wear latex or vinyl gloves while handling reagents and RNA samples to prevent RNase contaminations from surface of the skin or from dusty laboratory equipment.
- Change gloves frequently and keep tubes closed.
- Keep isolated RNA on ice.
- Reduce preparation time as much as possible.
- Use only sterile, disposable polypropylene tubes throughout the procedure (these tubes are generally RNase-free.)
- Non-disposable plastic ware should be treated before use to ensure that it is RNase-free. Plastic ware should be thoroughly rinsed with 0.1 M NaOH, 1 mM EDTA followed by RNase-free water. You can also take chloroform-resistant plastic ware rinsed with chloroform to inactivate RNases.

- All glassware should be treated before use to ensure that it is RNase-free. Glassware should be cleaned with detergent, thoroughly rinsed and oven baked at 240 °C for four hours or more before use. Autoclaving will not inactivate RNase activity completely. Oven baking inactivates RNases and ensures that no other nucleic acids (such as plasmid DNA) are present on the surface of the glassware. You can also clean glassware with 0.1 % DEPC (diethyl pyrocarbonate). The glassware has to be immersed in 0.1 % DEPC solution for 12 hours at 37 °C followed by autoclaving or heating to 100 °C for 15 minutes to remove residual DEPC.
- Electrophoresis tanks should be cleaned with detergent solution (e.g. 0.5 % SDS), thoroughly rinsed with RNase-free water, rinsed with ethanol and finally allowed to dry.
- All buffers have to be prepared with DEPC-treated RNase-free water.
- Avoid handling bacterial cultures, cell cultures or other biological sources of RNases in the same lab where the RNA purification will be performed.
- Do not use equipment, glassware and plastic ware employed for other applications which might introduce RNase contaminations in the RNA isolation.

4 Storage conditions

All kit components are shipped at ambient temperature.

Upon arrival store **MAG Suspension F** at 4 °C to 8 °C.

All other components of the kit should be stored dry at room temperature (15 °C to 30 °C). When stored at room temperature, the kit is stable until the expiration date printed on the label on the kit box.

If there are any precipitates within the provided solutions solve these precipitates by careful warming. Before every use make sure that all components have room temperature.

5 Functional testing and technical assistance

The IST Innuscreen GmbH guarantees the correct function of the kit for applications as described in the manual. This product has been produced and tested in an ISO 13485 certified facility.

We reserve the right to change or modify our products to enhance their performance and design. If you have any questions or problems regarding any aspects of the **innuPREP RNA Kit PP-Mini** or other IST Innuscreen GmbH products, please do not hesitate to contact us. For technical support or further information in Germany please contact info.innu@ist.com. For other countries please contact your local distributor.

6 Product use and warranty

The kit is not designed for the usage of other starting materials or other amounts of starting materials than those, referred to in the manual (→ "Product specifications" p. 9). Since the performance characteristics of IST Innuscreen GmbH kits have just been validated for the application described above, the user is responsible for the validation of the performance of IST Innuscreen GmbH kits using other protocols than those described below. IST Innuscreen GmbH kits may be used in clinical diagnostic laboratory systems after the laboratory has validated the complete diagnostic system as required by CLIA' 88 regulations in the U.S. or equivalents in other countries.

All products sold by IST Innuscreen GmbH are subjected to extensive quality control procedures and are warranted to perform as described when used correctly. Any problems should be reported immediately.

NOTE

This kit is for research use only!

7 Kit components

7.1 Components included in the kit

	Σ 16	Σ 96
REF	845-PS-0060016	845-PS-0060096
Lysis Solution RP	10 ml	50 ml
MAG Suspension F	0.25 ml	1.1 ml
Spin Filter	16	2 x 50
Receiver Tubes	20	2 x 50
Washing Solution HS (conc.)	7 ml	50 ml
Washing Solution LS (conc.)	6 ml	35 ml
RNase-free Water	5 ml	20 ml
Manual	1	1

7.2 Components not included in the kit

- 1.5 ml tubes
- 96 %–99.8 % ethanol (molecular biology grade, undenatured)
- Isopropanol (molecular biology grade)
- 96-well-plates or strips and tip combs for PurePrep Mini

7.3 Components needed for isolation from bacteria (not included)

- Lysozyme (stock solution 10 mg/ml (400 U/μl))
- Mutanolysin (stock solution 0.4 U/μl)
- Lysostaphin (stock solution 0.4 U/μl)
- 1 x TE-Buffer

Alternatively:

- innuPREP Bacteria Lysis Booster (845-KA-1000050, 50 rxn, IST Innuscreen GmbH)

7.4 Components needed for isolation from yeasts (not included)

- Yeast Digest Buffer (50 mM KH_2PO_4 , 10 mM DTT, pH 7.5)
- Lyticase (stock solution 10 U/ μl)

8 Product specifications

1. Starting material:

- Tissue sample (up to 20 mg)
- Eukaryotic cells (max. 5×10^6)
- Bacteria / Yeast cell pellets ($1 \times 10^8 - 1 \times 10^{12}$ cells)

2. Time for automated extraction protocol on PurePrep Mini:

- Approx. 20 minutes (excluding external lysis)

9 Initial steps before starting

- Add the indicated volume of absolute ethanol to **Washing Solution HS (conc.)** and mix thoroughly. Always keep the bottle firmly closed!

845-PS-0060016 Add 7 ml ethanol to 7 ml Washing Solution HS.

845-PS-0060096 Add 50 ml ethanol to 50 ml Washing Solution HS.

- Add the indicated volume of absolute ethanol to **Washing Solution LS (conc.)** and mix thoroughly. Always keep the bottle firmly closed!

845-PS-0060016 Add 24 ml ethanol to 6 ml Washing Solution LS.

845-PS-0060096 Add 140 ml ethanol to 35 ml Washing Solution LS.

- Avoid freezing and thawing of starting material.

10 Sample Preparation

10.1 Protocol 1: RNA extraction from tissue samples

NOTE

To maximize the final yield of total RNA a complete homogenization of tissue sample is important! For the homogenization of tissue samples, it is possible to use commercially available rotor-stator homogenizer. It is also possible to disrupt the starting material using mortar and pestle in liquid nitrogen and grind the tissue sample to a fine powder.

- A. **Homogenization of tissue samples using a rotor-stator homogenizer**
1. Transfer the weighed amount of fresh or frozen starting material in a suitable reaction vessel for the homogenizer.
 2. Add 450 µl **Lysis Solution RP**.
 3. Homogenize the sample.
 4. Transfer the homogenized tissue sample into a 1.5 ml reaction tube and place the sample in **Lysis Solution RP** for longer storage at -22 °C to -18 °C or use the sample immediately for isolation of total RNA following the protocol.

B. Disruption of the tissue sample using a mortar and pestle and liquid nitrogen

1. Transfer the weighed amount of fresh or frozen starting material under liquid nitrogen and grind the material to a fine tissue powder.
2. Transfer the powder into a 1.5 ml reaction tube. Don't allow the sample to thaw!
3. Add **450 µl Lysis Solution RP** and incubate the sample for appropriate time for a further lysis under continuous shaking for 30 minutes.
4. Finally place the sample under **Lysis Solution RP** for longer storage at -22 °C to -18 °C or use the sample immediately for isolation of total RNA following protocol.
5. After lysis spin down unlysed material by centrifugation at maximum speed for 1 minute.
6. Place a Spin Filter into a Receiver Tube. Transfer the supernatant of the lysed sample onto the Spin Filter. Centrifuge at 11,000 x g for 2 minutes. Discard the Spin Filter.

Do not discard the filtrate, because the filtrate contains the RNA!

7. Proceed with "Automated extraction using PurePrep Mini" on p.14.

10.2 Protocol 2: RNA extraction from eukaryotic cells

IMPORTANT

Please note that up to 5×10^6 cells can be processed.

1. Add **400 μ l Lysis Solution RP** to the cell pellet. Incubate for 2 minutes at room temperature. Resuspend the cell pellet completely by pipetting up and down. Incubate the sample for further 10 minutes at room temperature.
-

NOTE

To maximize the final yield of total RNA a complete disruption and lysis of the cell pellet is important! No cell clumps should be visible after lysis step.

4. Place a Spin Filter into a Receiver Tube. Transfer the lysed sample onto the Spin Filter D. Centrifuge at $11,000 \times g$ for 2 minutes. Discard the Spin Filter.

Do not discard the filtrate, because the filtrate contains the RNA!

5. Proceed with "Automated extraction using PurePrep Mini" on p.14.

10.3 Protocol 3: RNA extraction from bacteria and yeasts

IMPORTANT

We recommend a pre-incubation of bacterial cells with Lysozyme or optionally available **innuPREP Bacteria Lysis Booster**.

1. Collect the cells by centrifugation with parameters adequate for the cell type (e.g. 10 minutes at $3,000 \times g$). Discard the supernatant.
Do not discard the pellet!
2. Resuspend the bacterial cell pellet in **200 μ l TE Buffer** and the yeast cell pellet in **200 μ l Yeast Digest Buffer** (\rightarrow see "Components needed for isolation from yeasts", p. 9). in a 1.5 ml tube and add the

enzymes according to the table below or follow the instruction of the product innuPREP Bacteria Lysis Booster:

	Enzyme	Volume
Gram-negative bacteria	Lysozyme: 10 mg/ml (400 U/ μ l)	20 μ l
Staphylococcus spp.	Lysostaphin: 0.4 U/ μ l	10 μ l
Gram-positive bacteria	Lysozyme: 10 mg/ml (400 U/ μ l) + Mutanolysin: 0.4 U/ μ l	20 μ l + 5 μ l
Yeast	Lyticase: 10 U/ μ l	10 μ l

3. Incubate at 37 °C for 30 min under continuous shaking.
4. Add 200 μ l Lysis Solution RP to the pre-lysed sample and vortex vigorously or pipette sometimes up and down. Incubate the sample for further 10 minutes at room temperature.

NOTE

To maximize the final yield of total RNA a complete disruption and lysis of the cell pellet is important! No cell clumps should be visible after lysis step.

5. Place a Spin Filter into a Receiver Tube. Transfer the lysed sample onto the Spin Filter. Centrifuge at 11,000 x g for 2 minutes. Discard the Spin Filter.

Do not discard the filtrate, because the filtrate contains the RNA!

NOTE

If the solution has not completely passed through the Spin Filter, centrifuge again at higher speed or prolong the centrifugation time.

6. Proceed with "Automated extraction using PurePrep Mini" on p.14.

11 Automated extraction using PurePrep Mini

11.1 Prefilling of the DW Plate or the DW Strips

Cavity of KF96 DW Plate/Strip	Content
Cavity 1	400 µl lysed sample + 400 µl Isopropanol + 10 µl MAG Solution F
Cavity 2	800 µl Washing Solution HS
Cavity 3	800 µl Washing Solution LS
Cavity 4	800 µl Washing Solution LS
Cavity 5	empty
Cavity 6	100 µl – 150 µl RNase-free Water (volume depends on starting material and expected yield)

11.2 Loading filled Deep Well Plate/Strips to the PurePrep Mini and plug in the Tip Combs

NOTE

- When using strip (strips), the strip is inserted into the tray. In total, a maximum of 8 strips can be used in one extraction-run.
- The tip combs always dip staggered into the Strips.
Left tray side: Tip 1, 3,5, 7
Right tray side: Tip 2,4,6,8.
- It is recommended to mark the tips used for the extraction so that they are not used more than once

-
1. Select the protocol
"RNA1" and start the run.
 2. After finishing the extraction protocol, the Cavity 6 contains the isolated RNA.

3. Transfer the RNA into a fresh 1.5 ml Tube.

IMPORTANT NOTE

After finishing the extraction protocol, the last cavity of the Plate/Strip contains the isolated DNA/RNA. Store the DNA/RNA under adequate conditions.

We recommend storing the extracted RNA at $-82\text{ }^{\circ}\text{C}$ to $-78\text{ }^{\circ}\text{C}$.

If the eluate contains carryover of magnetic particles, place the plate on a magnet or centrifuge the plate at maximum speed for 3 minutes. Pipet the supernatant with DNA/RNA into a new plate.

12 Troubleshooting

Problem / probable cause	Comments and suggestions
Poor lysis of starting material	
Insufficient disruption or homogenization	After lysis centrifuge lysate to pellet debris and continue with the protocol using the supernatant. Reduce amount of starting material.
Little or no total RNA eluted	
Insufficient disruption or homogenization	Reduce amount of starting material. Overloading reduces yield!
DNA contamination	
Too much starting material	Reduce amount of starting material.
Incorrect lysis of starting material	Use the recommended techniques for lysis of cell pellet.
Total RNA degraded	
RNA source inappropriately handled or stored	Ensure that the starting material is fresh! Ensure that the protocol, especially the first steps, has been performed quickly.
RNase contamination of solutions; Receiver Tubes, etc.	Use sterile, RNase-free filter tips. Before every preparation clean up the pipette, the devices and the working place. Always wear gloves!

IST Innuscreen GmbH
Robert-Rössle-Str.10
13125 Berlin · Germany

Phone +49 30 9489 3380
Fax +49 30 9489 3381

info.innu@ist-ag.com